CONTEXTE:

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Total: pts/40 pts

SPECTR-O-MATIC QUICK REFERENCE GUIDE

1. GENERAL FEATURES OF THE SPECTR-O-MATIC:

The Spectr-o-Matic is a high-throughput spectrometer dedicated to routine analysis.

It allows automatic recording of absorbance spectra (Abs = 0-3 units) in the UV-visible region (200-800 nm), detection of unknown absorbing species (the library contains > 10'000 spectra), and determination of their concentration with a high level of accuracy.

The instrument is equipped with the following components:

- a set of UV and visible light sources (hydrogen lamp, respectively tungsten lamp)
- a quartz prism
- an automatically adjustable slit (0.2-2 nm)
- a compartment for a quartz cuvette containing the solution to be analysed
- a high-sensitivity photo-multiplier
- an electronic driver that controls the components
- a built-in computer that identifies unknown species and calculates concentrations

2. USE OF THE SPECTR-O-MATIC:

- a. Switch the power on (white button on the right side of the instrument); wait 10 min for warming up.
- b. Insert the quartz cuvette with the unknown solution in the cuvette compartment; close the cover lid.
- c. Start the automatic measurement (yellow button on the front panel); wait 30 sec for completion.
- d. Start the automatic identification (blue button on the front panel); wait 30 sec for completion.
- e. Check on the screen that the spectrum of the unknown species fits the spectrum of the identified species. In particular, check that the absorbances and the maximum wavelength (λ_{\max}) of the unknown and the identified species are the same.
- f. Print the results: spectrum, species name, species concentration (green button on the front panel).
- g. Repeat steps b-f for additional samples, if necessary.
- h. Remove the cuvette and switch the power off (black button on the right side of the instrument).

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
3. TROUBLESHOOTING:

In case of problems, do not use the repair kit! Call technical assistance (tel. 021-6933772, 24/24).

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Subject: order for spare parts

Miss, Mister,

I want to command the followed spare part for our instrument Spectr-o-Matic:

- 1 connexion cabel (part #101.17.23; \$18.35)
- 1 game of cuvettes for UV-absorbing solutions (part #101.52.19; \$55.90)
- 1 reparation kit: screwdriver, hamer, cork-screw (part #101.88.20; \$8.45)

I thank you four a quick delivery at the following adresse:

AnalyTeam S.A.

A. Tomik, lab technical

12, chemin des Rosiers

CH-1015 Lausanne

Best regards.